


BioShipper-2, for UN3373 Biological Substances

Conforms with Packaging instruction 650




BioBag 752
230 x 315mm
Int Dims
(Code 752)

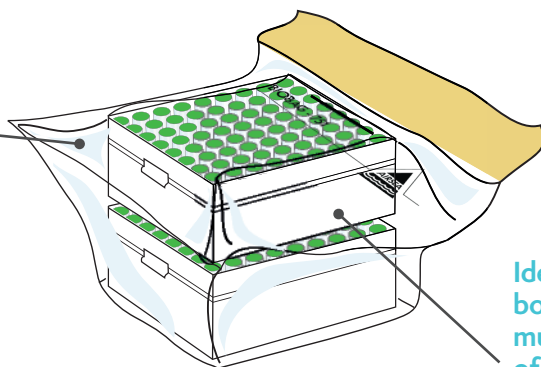

Ideal for cryovial
boxes or a
multitude
of single primary
containers
(Not Supplied)

BioShipper-2
315 x 273 x 101mm
Int Dims
(Code 762)

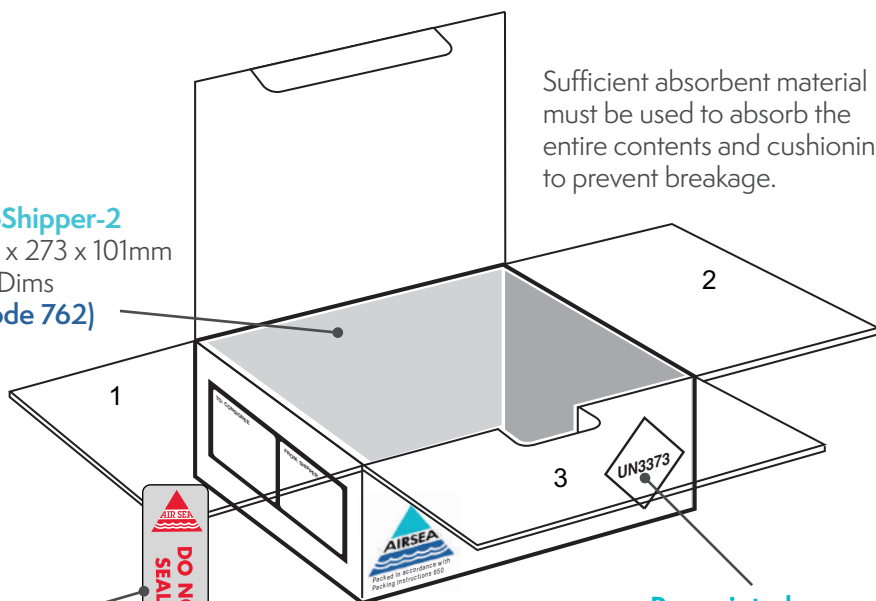

**Foil Tamper
Evident Seal
(Code 741)**



**Pre-printed
UN3373 Label**

Add additional cushioning to fill void spaces

ASSEMBLY DETAILS: **BioShipper-2** Code: 762

Box Dimensions:
External 318 x 276 x 110mm
Internal 315 x 273 x 101mm

Instructions:

Place primary receptacle(s) into the cushioning pouch(s). If using cryovial use absorbent sheet

Place into the 95kPa bag and seal as per printed instructions.

Place sealed bag(s) into BioShipper-2. Seal with tamper evident seal.

Notes:

The BioShipper-2 is supplied flat packed, it's designed with an easy pop up construction. Capable of carrying up to 2 cryovial boxes in 1x BioBag 752, or multiple single primary receptacles in numerous biobag 752.

The BioShipper-2 comes complete with:

Code 752, 95Kpa BioBag

Code 743, Security Seal

Code 740, 4 Bay Absorbent
Cushioning Pouch x 2 (240ml)

Outer Box with Pre-Printed
UN3373 Label

Thermal control:

A variety of Thermal Control Units for transportation in dry ice or coolpack are available.

It is the Shippers responsibility to ensure that packagings conform to the relevant transport regulations.

Further info



CAUTION

This packaging is NOT a U.N. Certificated Packaging for the transport of class 6.2 Infectious Substances, Category A and does not conform to Packaging Instructions 620.

Air Sea Containers Ltd

521 Cavendish Street, Merseyside, CH41 8FZ, UK

▲ Tel +44 (0)151 653 1500 ▲ Email uk@airseadg.com ▲ Visit www.airseadg.com